TR

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EN

The newly developed jaw crusher rotary sample divider is an excellent choice for industrial and laboratory uses. The product takes materials with an inlet size of < 100 mm and manages to crush them down to an exit fineness of 0-20 mm. This superior crushing capacity can process up to 150 kg of material per hour, making it ideal for high volume testing needs. The sample splitting feature is distinguished by the ability to separate $\pm 1\%$ into 8 identical representative samples by weight. This feature increases the accuracy and reliability of test results, so more precise analysis can be made.

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EN

The device has 8 4-litre compartments and can separate samples into eight equal representative samples of $\pm 1\%$ by weight. This feature is especially essential for comprehensive analyzes and quality control processes. Precision and consistency were prioritized during the sample division process.

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ROTARY SAMPLE DIVIDER WITH JAW CRUSHER**About This Performance Data**

This sheet shows how well our machines work. You'll see numbers for how much they can handle, how fast they work, and the size of the stuff they make. It's all about helping you use our machines the best way.

Rotary Sample Divider with Jaw Crusher	
Sample & Size	Marble 20 mm
Sample Weight	123 kg
Gap between Jaws (output)	3 mm
Divider Speed	25 -30 rpm
Crushing & Dividing Time	35 minutes
Capacity	210 kg/h

DIVIDER DISTRIBUTION			
Divisions	Net Weight (kg)	Mean Difference (kg)	Mean Difference (%)
Division 1	15,30	-0,1	-0,4
Division 2	15,22	-0,1	-0,9
Division 3	15,46	0,1	0,7
Division 4	15,22	-0,1	-0,9
Division 5	15,49	0,1	0,9
Division 6	15,51	0,2	1,0
Division 7	15,44	0,1	0,5
Division 8	15,23	-0,1	-0,8
Total Weight	122,87		
Mean	15		